
Product Specification

Product Name: FB2 Additives
Product Code: SG7291 / SG7292 / FB2-set

PRODUCT DESCRIPTION:

Acts as a supplement for FB2 (Full-Fraser).

USE:

Before opening the ampoule, disinfect the surface with a 75% alcohol swab. Open the ampoule under sterile conditions. Add one vial each of FB2 Additive A and B or the FB2-set (FB2 Additives) to every 100 mL of Fraser Broth Medium.

Item	Specification
Packing Size:	1 mL x 10 Vials/box
Formulation (Reagent A):	2.0 mg Naphthylpyridinone Acid, 2.5 mg Acriflavin hydrochloride
Formulation (Reagent B):	0.05 g Ferric ammonium citrate
Storage Condition:	2-8°C
Shelf Life:	1 Year

Waste Disposal: After usage, bacteria-contaminated items should be autoclaved at 121°C for 30 minutes.